

**Advance the potential of your cell culture**  
*Cedex Bio and Cedex Bio HT Analyzers*



## The value of certainty

### *Cedex Bio and Cedex Bio HT Analyzers*

#### *Understanding the challenges of bioprocess monitoring*

Whether you are working in the pharmaceutical or biotech industry, you always value the safety and confidence of monitoring and analyzing cell culture with state of the art solutions.


As one of the leaders in the diagnostics and pharmaceutical field, Roche has first-hand experience with challenges in biomanufacturing. That knowledge, plus many years of competence in designing diagnostic tests and engineering analyzer systems, are the foundation for the Cedex Bio Analyzer family from Roche CustomBiotech.

The Cedex metabolite analyzers are based on enzymatic assays, enabling high robustness and sensitivity as well as optimized measurement ranges, adapted to the particular demands of cell culture and fermentation processes. Thus, you can rely on one instrument family to capture useful and timely insights into your bioprocess and make the right decisions for optimal manufacturing outcomes.

We continuously improve the performance and level of integration of our analyzers to support you with reliable and up-to-date systems. Moreover we offer a comprehensive test menu which we are constantly evolving according to the user needs, so you can flexibly monitor multiple and modern parameters in your bioprocess without having to develop new assays on your own or purchase additional instruments.

#### *Act on predictability and accuracy*

Ultimately, having a bioprocess under control means predicting changes in cell cultivation and fermentation that can impact product yield and quality. This is especially true when running a bioprocess within narrow specifications, such as under substrate-limiting conditions. High predictability requires precise and accurate data that reflect the actual conditions in a bioreactor, even at low analyte concentrations. Compared to membrane-based analyzers, the Cedex Bio Analyzers offer significantly higher reproducibility, accuracy and precision (Figure 1). For example, you can rely on the monitoring results to indicate the right time points for substrate bolus shots or harvest.

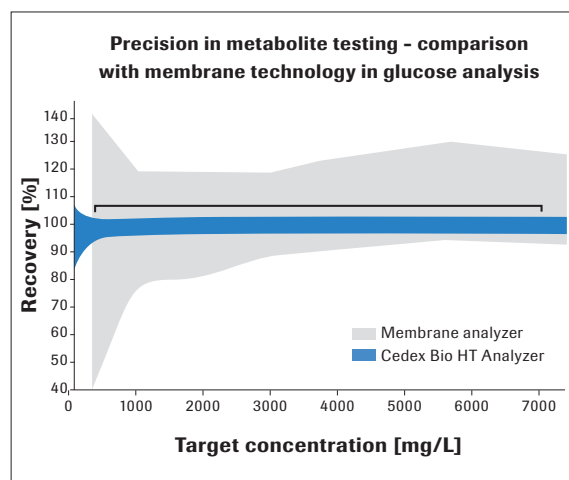


**Figure 1:** Recovery of glucose concentration was determined in parallel using the Cedex Bio HT Analyzer and a membrane analyzer. (Roche evaluation data)

### *See relevant information in time*

Information about the conditions in a cell culture must be available in time to enable controlled processing with optimal manufacturing outcome. Thus, the Cedex Bio Analyzers offer an assortment of assays, some designed specifically to determine the product yield and quality throughout the fermentation process. Quantify IgG levels in minutes and with the same accuracy as the gold standard HPLC (Figure 2).

In addition, determine lactate dehydrogenase (LDH) as a sensitive control for cell integrity. A rising LDH activity in the culture medium is an early indicator of cell lysis, which is combined with the release of intracellular proteases, glucosidases and other degrading proteins, generating a considerable risk for the yield and quality of the valuable product.

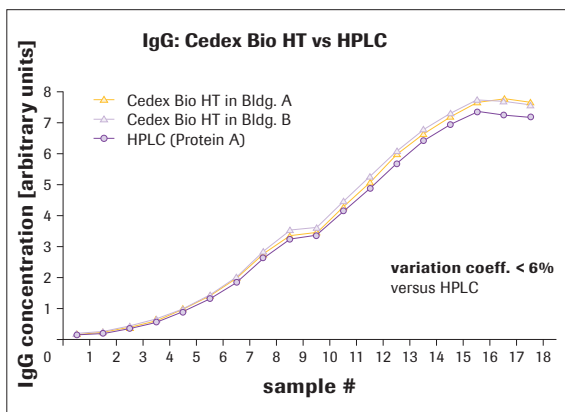

**Figure 2: Results comparable to HPLC.** The IgG titer in the culture medium was monitored over a production period of 18 days. Samples were analyzed in parallel using HPLC (Protein A) and two Cedex Bio HT Analyzers in separate locations (bldg A/B). The consistent results represent the perfect precision and the equivalence of the methods. (Roche evaluation data)

### *Count on answers, every time*

Cedex Bio Analyzers were developed based on industry-proven diagnostic technology. Adapted to support process development and large-scale production at Roche Pharmaceuticals, each instrument withstands the demanding conditions in the daily bioprocessing, with need for high robustness and minimal manual interaction.

**The flexible capabilities for automated sample pre-dilution, the long on-board stabilities of the reagents and the long calibration intervals, the dedicated multi-analyte calibrators and controls, and the identification of reagents and samples via barcodes, make Cedex Bio Analyzers a perfect match for highly efficient process monitoring.**

## Powerful, flexible and comprehensive cell culture profiling

### *Cedex Bio and Cedex Bio HT Assays*

*Find answers you need, with a comprehensive testing panel*

Each available assay facilitates tight control of biopharmaceutical manufacturing, and new assays are consecutively added to the menu based on industry knowledge and requests from customers. Thus, with a single platform, you have continuously increasing testing flexibility and insights to optimize your bioprocess.



Additional assays are in development.

Visit [custombiotech.roche.com/cedex](http://custombiotech.roche.com/cedex) for updates or contact your local CustomBiotech team.

# Select the right analyzer to optimize your bioprocess workflow

*Different specifications, identical test results*




Cedex Bio Analyzer

Cedex Bio HT Analyzer

Throughput	
60 – 80 tests per hour	260 – 400 tests per hour
Sample capacity	
up to 8 samples at once (continuous loading)	up to 96 + 30 samples at once (continuous loading)
Test kit capacity	
13	32
Daily maintenance	
semi automated, <15 – 20 minutes hands on	automated
Automated on-board sample dilution	
selectable dilution	freely programmable

*Cedex Bio and Cedex Bio HT Analyzers deliver consistent results*

Although the two analyzers differ in throughput and certain aspects you can rely on equivalent results. High correlation between the Cedex Bio and the Cedex Bio HT Analyzers has been proven for all assays.

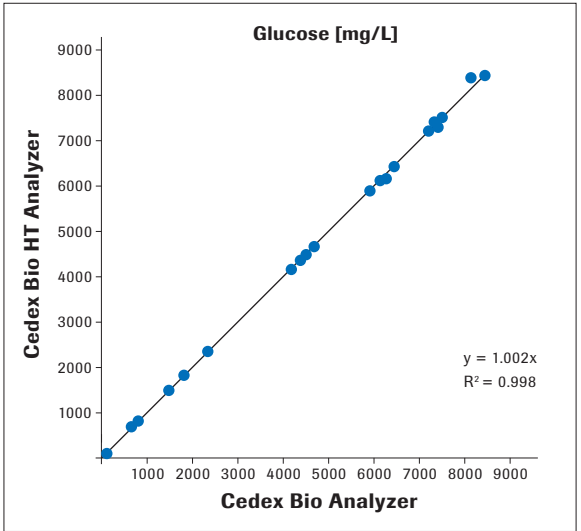


**Figure 3:** Comparison of the Cedex Bio and Cedex Bio HT Analyzer showing the same recovery of glucose. (Roche evaluation data)



## Cedex Bio Analyzer

*Small footprint, big impact*

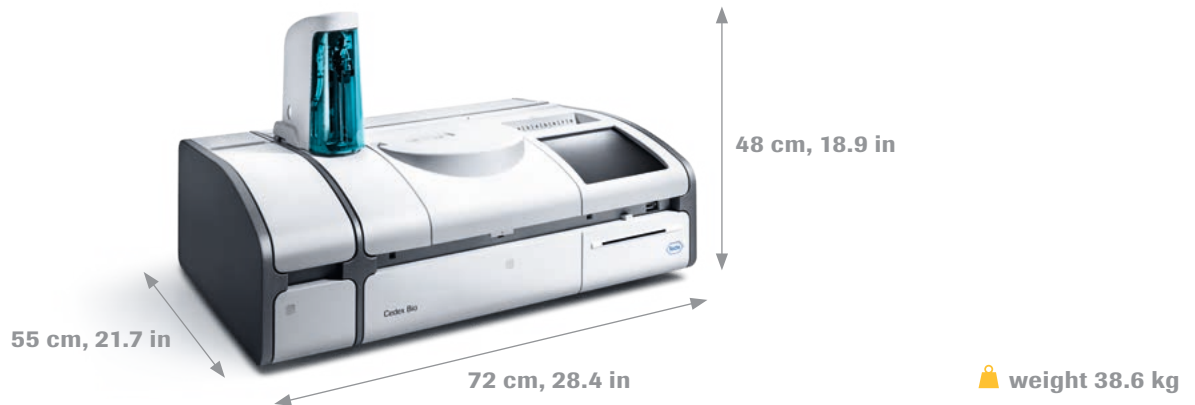

*A reliable, high-quality instrument for small sample throughput*

- **Small size:** Benchtop system suitable for any space
- **Easy handling:** Sophisticated and intuitive operating software
- **Integrated processing unit:** Built-in touchscreen reduces qualification effort
- **Reliable answers:** Robust and proven technology
- **Efficiency:** Low maintenance effort and frequency save time and money

### Cedex Bio and Bio HT Analyzer

- Same menu of assays with high specificity and sensitivity
- LIMS connectivity through middleware systems with complete traceability and regulatory compliance


## Cedex Bio HT Analyzer

*High efficiency, fast action*



*The right instrument for improved workflow efficiency*

- **High throughput:** Numerous samples can be processed simultaneously with minimal manual intervention
- **Enhanced flexibility with the 96-well Upgrade Kit:**
  - Use 96-well plates or conventional racks (total capacity: up to 96 + 30 samples)
  - Boosts processing safety and efficiency, minimizing overall hands-on time (e.g., 96 glucose tests in 35 min)
  - Designed for manual or robotic loading of 96-well plates
- **Long-term operation:** Long time on-board stability of reagents and long assay calibration intervals
- **Flexible dilution:** Automated sample dilution capabilities save time and ensure high accuracy
- **Full automation:** Online integration with automated systems to optimize measurement points and timelines

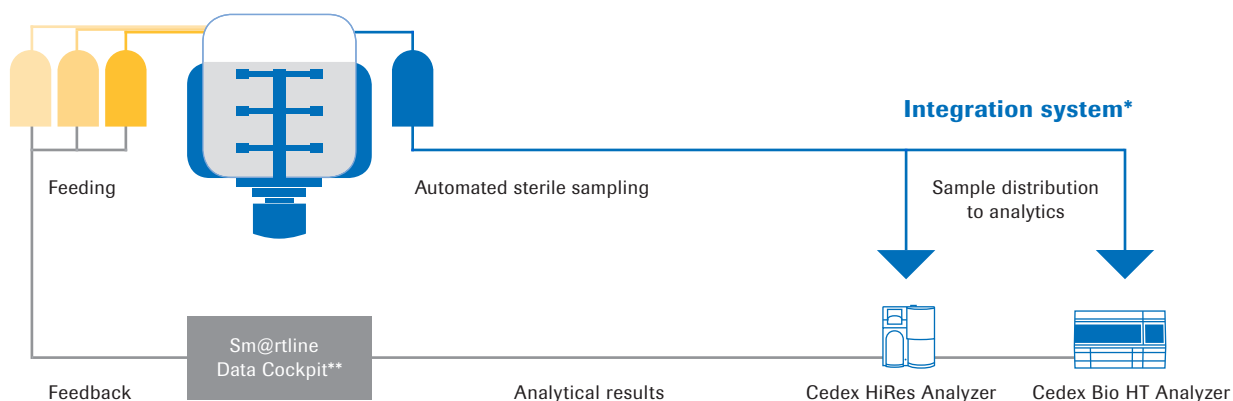

**Figure 4:** Process integration and automation capabilities with Cedex Bio HT and Cedex HiRes Analyzer

\*Various integration systems are available.

\*\*Sm@rtline Data Cockpit is provided by AGU, Leverkusen, Germany

## Ordering information

Product	Catalog number
Cedex Bio Analyzer	06 395 554 001
Cedex Bio HT Analyzer	06 608 116 001
Kit 96-well plate Cedex	09 181 687 001
Rack 96-well plate Cedex	09 241 400 001

The ordering information for all Cedex Bio Analyzer and Cedex Bio HT Analyzer assays is available on **[custombiotech.roche.com](https://custombiotech.roche.com)**

## Regulatory disclaimer

For use in quality control/manufacturing process only.

## Trademarks

CEDEX is a trademark of Roche.

**[custombiotech.roche.com/cedex](https://custombiotech.roche.com/cedex)**

## Please contact your local CustomBiotech representative

### Europe, Middle East, Africa, Latin America

Phone +49 621 759 8580  
Fax +49 621 759 6385  
[mannheim.custombiotech@roche.com](mailto:mannheim.custombiotech@roche.com)

### United States

Phone +1 800 428 5433, ext. 14649 (toll-free)  
Fax +1 317 521 4065  
[custombiotech.ussales@roche.com](mailto:custombiotech.ussales@roche.com)

### Canada

Phone +1 450 686 7050  
Fax +1 450 686 7012  
[custombiotech.can@roche.com](mailto:custombiotech.can@roche.com)

### Japan

Phone +81 3 6634 1046  
Fax +81 3 5479 0585  
[japan.custombiotech@roche.com](mailto:japan.custombiotech@roche.com)

### Asia Pacific

Phone +65 6371 6638  
Fax +65 6371 6601  
[apac.custombiotech@roche.com](mailto:apac.custombiotech@roche.com)

### Published by

Roche Diagnostics GmbH  
Sandhofer Straße 116  
68305 Mannheim  
Germany

© 2020 Roche Diagnostics GmbH  
All rights reserved.